


Rapid immunological detection systems

pab ELISA-Stick with test reagent set


Pab ELISA-Stick was designed for qualitative analysis of antigenic substances in soluble and insoluble samples. The pab ELISA-Stick Assay is a sensitive and easy to handle test for the detection of various substances. The test can be carried out without sample preparation or pre-incubation. Since the test does not need any apparatus equipment, it is especially suitable for analysis on the spot.

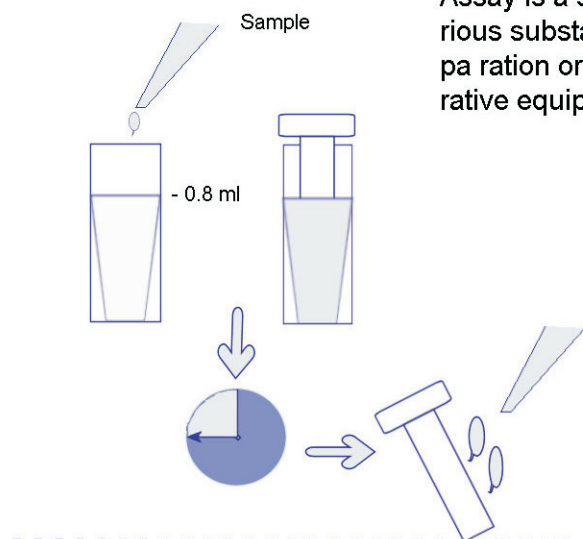

Fill vial 2 with the sample up to 0.8 ml.
Stick the pab ELISA-Stick in vial with the sample.

Incubate 45 min at room temperature.


Wash the ELISA-Stick with water.

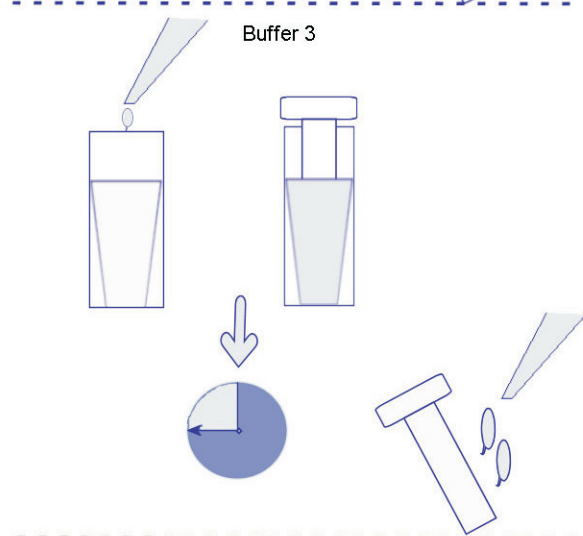

Fill vial 4 with buffer from vial 3.
Stick the pab ELISA-Stick in vial 4.

Incubate 45 min at room temperature.

Wash the ELISA-Stick with water.

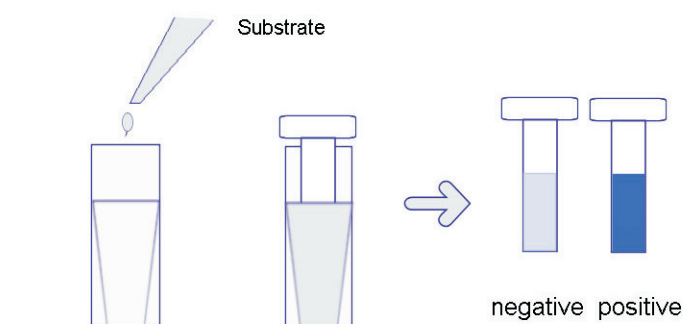

Fill the substrate solution from vial 5 into the cuvette.
Stick the pab ELISA-Stick into the cuvette.

Expect the colour reaction.

Rapid immunological detection systems

Product name	Product No.	Sensitivity	Package size
anti-aflatoxin M1 ELISA-Stick	EA-2	1 ng	5 Tests
anti-aflatoxin M1 ELISA-Stick	EA-3	1 ng	10 Tests
anti-aflatoxin B1 ELISA-Stick	EA-4	1 ng	5 Tests
anti-aflatoxin B1 ELISA-Stick	EA-5	1 ng	10 Tests
anti-Aspergillus niger ELISA-Stick	EA-6	1 ng	5 Tests
anti-Aspergillus niger ELISA-Stick	EA-7	1 ng	10 Tests
anti-chloramphenicol ELISA-Stick	EC-2	1 ng	5 Tests
anti-chloramphenicol ELISA-Stick	EC-3	1 ng	10 Tests
anti-chlorinated phenols ELISA-Stick	EP-8	1 ng	5 Tests
anti-chlorinated phenols ELISA-Stick	EP-9	1 ng	10 Tests
anti-deoxynivalenol ELISA-Stick	ED-2	1 ng	5 Tests
anti-deoxynivalenol ELISA-Stick	ED-3	1 ng	10 Tests
anti-dinitro-aromates ELISA-Stick	ED-5	1 ng	5 Tests
anti-dinitro-aromates ELISA-Stick	ED-6	1 ng	10 Tests
anti-ergotamine ELISA-Stick	EE-2	1 ng	5 Tests
anti-ergotamine ELISA-Stick	EE-3	1 ng	10 Tests
anti-Gaeumannomyces graminis ELISA-Stick	EG-2	1 ng	5 Tests
anti-Gaeumannomyces graminis ELISA-Stick	EG-3	1 ng	10 Tests
anti-ochratoxin A ELISA-Stick	EO-2	1 ng	5 Tests
anti-ochratoxin A ELISA-Stick	EO-3	1 ng	10 Tests
anti-patulin ELISA-Stick	EP-6	1 ng	5 Tests
anti-patulin ELISA-Stick	EP-7	1 ng	10 Tests
anti-penicillin G ELISA-Stick	EP-1	1 ng	5 Tests
anti-penicillin G ELISA-Stick	EP-2	1 ng	10 Tests
anti-penicillin B ELISA-Stick	EP-4	1 ng	5 Tests
anti-penicillin B ELISA-Stick	EP-5	1 ng	10 Tests
anti-phenyl urea ELISA-Stick	EP-10	1 ng	5 Tests
anti-phenyl urea ELISA-Stick	EP-11	1 ng	10 Tests
anti-Ramularia ELISA-Stick	ER-6	1 ng	5 Tests
anti-Ramularia ELISA-Stick	ER-7	1 ng	10 Tests
anti-rubellin ELISA-Stick	ER-4	1 ng	5 Tests
anti-rubellin ELISA-Stick	ER-5	1 ng	10 Tests
anti-saxitoxin ELISA-Stick	ES-8	1 ng	5 Tests
anti-saxitoxin ELISA-Stick	ES-9	1 ng	10 Tests
anti-triazine ELISA-Stick	ET-5	1 ng	5 Tests
anti-triazine ELISA-Stick	ET-6	1 ng	10 Tests
anti-tylosin ELISA-Stick	ET-2	1 ng	5 Tests
anti-tylosin ELISA-Stick	ET-3	1 ng	10 Tests
anti-zearalenone ELISA-Stick	EZ-1	1 ng	5 Tests
anti-zearalenone ELISA-Stick	EZ-2	1 ng	10 Tests